

TITLE: SuperScript III First Strand cDNA synthesis

SOP Number: D-DEZ-PRO-019

Revision Number: 0

Effective Date: 07 Apr 2015

6. Terminate reactions at 85 degrees Celsius for 5 min. Chill on ice.
7. Add 1 µl RNase H to each sample and incubate at 37 degrees Celsius for 20 min.
8. cDNA samples can be stored at -20 degrees Celsius or used immediately for PCR.

F. HISTORY

Effective Date	Revision	Change
07 Apr 2015	0	Initial issue of SOP