

TITLE: Logging onto Iowa Courses ON-line (ICON) for GMP Related Training

SOP Number: D-DEZ-EQP-001

Revision Number: 0

Effective Date: 07 Apr 2015

and will contact Quality Assurance to order an additional Dewar. (The turn-around can be as soon as 24 hours)

1.2 Troubleshooting

1.2.1 Power Supply

1. The ProStar panel LED indicator is not sensing an alarm when the Dewar is empty:
 - Check that the cord is securely plugged in.
 - Plug another appliance into the outlet to see if it is live.
 - If the outlet is dead, check the circuit breaker or fuses.
 - Check the Pressure Switches are operating correctly.

1.3 Alarms

- 1.3.1 The Prostar Alarm system is designed to provide a signal when the Nitrogen Dewar is empty. The alarm will send notification to the Johnson Controls MVE and to the system owner.
- 1.3.2 The alarm is activated when the “Indicating Pressure Switch” detects a pressure loss and works in conjunction with the ProStar alarm panel. The audible warning signal sounds when there is a pressure loss condition.
- 1.3.3 Standard operating conditions of the two-point alarm are indicated by an illuminated green lamp. The lamp and horn check button allows for self-checking.
- 1.3.4 If one or more alarm signals are triggered, acoustic and visual signals are emitted.

E. HISTORY

Effective Date	Revision	Change
07 Apr 2015	0	Original document