



# The University of Iowa Dezii Translational Vision Research Group

**TITLE: Preparation of Mounting Media**

SOP Number:           D-DEZ -PRO-010          

Revision Number:           0          

Effective Date:           07 Apr 2015          

---

5. If desired add DAPI at 2 ug/ml.
6. Aliquot the supernatant (we use 1 ml aliquots which is enough for 15-20 slides) and store at -20°C.
7. Since it polymerizes upon contact with air you should not refreeze the aliquots. Thaw it just before use and throw away what you do not use.

G. HISTORY

Effective Date	Revision	Change
07 Apr 2015	0	PM-18 Rev. #3